

# BioProcess International

BIOPROCESS TECHNICAL


## An Approach to Generating Better Biosimilars

### Considerations in Controlling Glycosylation Variability in Protein Therapeutics

Lori Fortin, Jonathan Fura, Hong Li, Brian Beyer, and Nandu Deorkar

The global market for biotherapeutics has expanded extensively over the past decade and is projected to account for more than a quarter of the pharmaceutical market by 2020, with sales exceeding US\$290 billion (1). Continued expansion of the biosimilar marketplace has led to many commercial opportunities and technical challenges. The biological systems used to manufacture such drug products are inherently variable

**Figure 1:** Different protein glycoforms; recombinant proteins expressed for use as biosimilars might not be FDA approved because of variation in protein glycosylation patterns that potentially alter protein structure, stability, and activity.

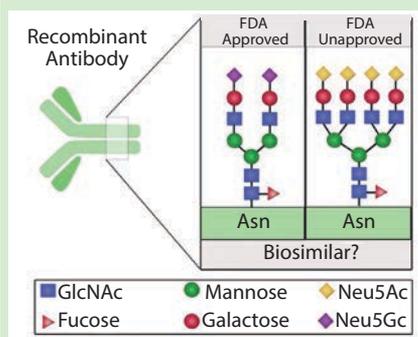

— a feature that has important consequences for the reproducibility, safety, and efficacy of the resulting products. Therefore, a prerequisite for introducing such biologics into routine clinical use is to ensure consistency of lot quality. Process understanding and consistency are critical because slight changes can lead to adverse effects

such as immunogenicity, with potentially serious safety implications. In generating biosimilars, it is crucial to ensure that all clinically active attributes of the biotherapeutic, including protein posttranslational modifications (PTMs), are accounted for and confirmed (2).

One such PTM that must be carefully monitored is protein glycosylation, which is the attachment of carbohydrates to a protein structure. That occurs within either the endoplasmic reticulum (ER) or the Golgi apparatus of a cell (GA) (3–5). Because those glycosylation mechanisms depend on factors such as cellular host, glycotransferase expression, substrate specificity, and substrate availability, different glycoforms of a single protein can be generated (3). Protein glycosylation can be a critical quality attribute for proteins used in discovery and biotherapeutics because it often plays a key role in their efficacy, safety,

**PRODUCT FOCUS:** PROTEINS

**PROCESS FOCUS:** PRODUCTION

**WHO SHOULD READ:** PROCESS DEVELOPMENT

**KEYWORDS:** GYCOFORMS, RECOMBINANT ANTIBODIES, BIOEQUIVALENCE, CULTURE MEDIA

**LEVEL:** INTERMEDIATE

stability, and functionality (6–10). Thus, regulatory agencies stress limited variability and note the importance of maintaining uniform glycosylation patterns during manufacturing and scale-up as well as for bioequivalence studies. As the market for protein biologics and biosimilars continues to expand, the need for generating homogeneous protein glycoforms through recombinant technology is of paramount importance (Figure 1). Herein, we identify possible strategies to mitigate glycosylation variability that can be used in generation of uniform proteins for discovery, therapeutics, and bioequivalents.

### CELL LINE VARIATION

During the discovery phase of a process, one of the first considerations for biotherapeutic protein production is the choice of host cell. Several host systems can be used, including *Escherichia coli*, yeast, and insect and mammalian cells. Ultimately, the choice depends on the complexity of a protein and required structural components to ensure activity, stability, efficacy, and quality. A number of review papers cover the individual benefits of each host system (11–14).

During discovery, several hosts can be screened to determine which will offer better protein yield, functionality, recoverability and overall quality. Organisms such as *E. coli* express proteins that have relatively simplistic folding requirements (e.g., lack of disulfide bridge formation, carbohydrate attachment, and lack of cofactor processing events). Although *E. coli* serves a limited role in the common production of proteins for discovery and commercial development, this system provides the unique opportunity of overexpressing quantities of raw polypeptide that can be refolded with varying results.

As an alternative, generation of biologically active proteins that require in vivo folding can be produced in yeast and baculovirus-infected insect cells (15). Yeast is capable of PTMs but typically results in high-mannose N-glycosylation (in the form of hyperglycosylation) and not complex-

ordered glycan attachment. If the glycosylation profile is not critical, yeast often can be a cost-effective and faster-growing and faster-producing host. However, yeast produces high levels of proteases, which can adversely affect product yield through protein degradation (16).

Plant systems are being considered for simple proteins and those produced using glycoengineering (12, 17). Plants can produce proteins in a short time, but they cannot perform complex glycosylation unless glycoengineering is used to express proteins similar to those produced in humans.

In attempts to generate proteins for discovery research and biotherapeutic application, mammalian cell recombinant expression can produce proteins that correctly mimic properties found within a human patient. Traditionally, mammalian cells have been the most commonly used hosts for producing glycosylated proteins such as monoclonal antibodies (MAbs) and enzymes, with many licensed products expressed by Chinese hamster ovary (CHO) cells. Because mammalian cells are more likely to produce proteins that are properly folded and contain the correct PTMs, the use of such cells presents several advantages over the alternative hosts discussed above. However, production from mammalian cells often is hindered by slower growth rates, lower protein yields, and higher costs of production. In addition to CHO, other mammalian cells used often include human embryonic kidney (HEK 293), myelomas such as NS0 mouse myeloma, hybridomas, baby hamster kidney (BHK), and PER. C6 human retinal cells (15). Because the origins of these cell lines vary, they can exhibit differences in cellular growth, protein production levels, and glycosylation patterns that must be considered by companies selecting host cells.

Further complicating the choice of host system, cell lines generated from clones can exhibit different glycosylation patterns. Therefore, several clones will be generated during discovery using a chosen host(s) and screened for protein production and quality (14, 18–22). As the fields of biosimilars and therapeutic proteins

grow, regulatory agencies are placing greater emphasis on controlling critical protein attributes such as glycosylation patterns. Therefore, when choosing a clone, it is necessary to ensure that the correct glycosylated protein is produced, even if production levels are not optimum. It is also beneficial to confirm that production in constitutive expression remains stable over the course of many generations, not only while establishing cell lines from individual clones, but also subsequently during development of the final manufacturing process.

### MEDIA SUPPLEMENTATION

Components within a medium can affect protein PTMs (9). One method that has been investigated to gain greater control of protein glycosylation is supplementation of the media with diverse small molecules such as sugars, amino acids, and trace elements. For example, media supplementation with 25 mM glucose (a main energy source for cells) enabled production of chimeric MAbs exhibiting increased rates of glycosylation and greater homogeneity of the protein glycoforms (23).

Different studies have been performed to regulate growth, protein production, and accumulation of lactate in cell cultures by using alternative sugars such as galactose and fructose, which are slowly metabolized. Using those sugars also can affect protein glycosylation. In one study, researchers found that complete substitution of glucose with galactose or fructose resulted in high-mannose forms of glycan in CHO cultures producing IgG (24). As expected, when cells were fed glucose and galactose simultaneously, that percentage dropped. In a second study, supplementation of media with monosaccharides (such as galactose and glucosamine) and the nucleotide uridine led to drastic increases in the sialylation of human cytokine interferon gamma when recombinantly expressed in CHO cells (25).

Similar to augmenting the sugar concentration within media, spiking amino acids into the media can affect recombinant protein glycosylation

greatly (26). Cellular respiration leads to the accumulation of ammonia, which can be toxic to cells and can alter recombinant protein glycosylation (27–29). To reduce ammonia generation, researchers often substitute glutamine with glutamate (30). And supplementation of media with proline, threonine, and glycine can mitigate the toxic effects of ammonia production and result in increased sialic acid and *N*-acetylglucosamine content of recombinantly expressed tissue plasminogen activator (t-PA), a protein drug used to treat stroke patients (29, 31). In a separate study, addition of asparagine to cell culture media altered the glycoforms of recombinant proteins. Such glycoform variation has been hypothesized to result from the need for asparagine in N-linked glycosylation and mitigation of toxic ammonium effects within cellular metabolism (32).

Addition of cofactors such as metal ions also can affect protein glycosylation significantly. Metals such as iron, cobalt, and copper have been added to cell culture media and were all shown to alter glycosylation levels of diverse proteins in unique cell lines (33–35). Similarly, supplementation of culture media with exogenous manganese can increase sialylation of human erythropoietin, a protein hormone used to treat anemia. That escalation in protein sialylation increased protein half-life within human serum and enhanced efficacy (36). Furthermore, the addition of manganese chloride along with uridine and galactose to cell culture media drastically increased the galactosylation of recombinant antibodies produced in CHO cells. Those results indicate that supplementation of cell culture media with certain supplements can have synergistic effects in enhancing production of a particular protein glycoform (37). Using that information, researchers demonstrated that galactosylation of recombinant monoclonal antibodies expressed in unique CHO cell lines can be controlled using those additives in specific ratios (38).

Further amplifying the effects of

cofactors present within a medium, trace concentrations of metals can potentiate protein glycoforms. One study determined that manganese impurities from different ferrous sulfate lots generated different glycosylation profiles of MAbs (39). In a separate study investigating the effects of iron and manganese on protein glycosylation, researchers observed that parts-per-million (ppm) concentrations of iron were required for altering protein glycosylation, whereas manganese altered protein glycoforms at concentrations of 20 parts per billion (ppb) (40). Such studies highlight the importance of using highly defined materials when producing media formulations because even trace impurities can produce undesirable effects on protein glycosylation.

#### KEY TO BIOSIMILARS APPROVALS

As the use of biotherapeutics continues to expand within the pharmaceutical marketplace, the development of biosimilars represents great potential for vast expansion within the industry. Imperative to the use and approval of these biosimilars is the understanding of protein glycosylation and the production of particular glycoforms. Although the generation of specific protein glycoforms remains poorly understood, the limited scientific data that we have indicate the need for a highly customized approach when producing biosimilars through recombinant technology. Production of precise protein glycoforms will demand specific cell lines and unique cell culture media formulations consisting of exceptionally defined reagents. With the market of biosimilars greatly expanding within the next decade, understanding protein glycosylation variability and the requirements necessary for generating specific biosimilars will be imperative to gain FDA approval and to address the needs of this new era of pharmaceuticals.

#### REFERENCES

1 Jacoby R, et al. *Winning with Biosimilars: Opportunities in Global Markets*.

Deloitte Development LLC: New York, NY, 2015.

2 Bolton A. AbbVie's Open Public Hearing Comments at FDA Arthritis Advisory Committee Meeting, 9 February 2016.

3 Kornfeld R, Kornfeld S. Assembly of Asparagine-Linked Oligosaccharides. *Annu. Rev. Biochem.* 54, 1985: 631–664; doi:10.1146/annurev.bi.54.070185.003215.

4 Van den Steen P, et al. Concepts and Principles of O-Linked Glycosylation. *Crit. Rev. Biochem. Mol. Biol.* 33(3) 1998: 151–208.

5 Breitling J, Aebi M. N-Linked Protein Glycosylation in the Endoplasmic Reticulum. *Cold Spring Harb. Perspect. Biol.* 5(8) 2013: a013359; doi:10.1101/cshperspect.a013359.

6 Abes R, Teillaud JL. Impact of Glycosylation on Effector Functions of Therapeutic IgG. *Pharmaceuticals* (Basel) 3(1) 2010: 146–157.

7 Sola RJ, Griebenow K. Effects of Glycosylation on the Stability of Protein Pharmaceuticals. *J. Pharm. Sci.* 98(4) 2009: 1223–1245; doi:10.1002/jps.21504.

8 Wright A, Morrison SL. Effect of Glycosylation on Antibody Function: Implications for Genetic Engineering. *Trends Biotechnol.* 15(1) 1997: 26–32.

9 Bruhlmann D, et al. Tailoring Recombinant Protein Quality By Rationale Media Design. *Biotechnol. Progress* 31(3) 2015: 615–629; doi:10.1002/btpr.2089.

10 Sha S, et al. N-Glycosylation Design and Control of Therapeutic Monoclonal Antibodies. *Trends Biotechnol.* 34(10) 2016: 835–846.

11 Butler M, Spearmann M. The Choice of Mammalian Cell Host and Possibilities for Glycosylation Engineering. *Curr. Opin. Biotechnol.* 30, 2014: 107–112; doi:10.1016/j.copbio.2014.06.010.

12 Durocher Y, Butler M. Expression Systems for Therapeutic Glycoprotein Production. *Curr. Opin. Biotechnol.* 20(6) 2009: 700–707; doi:10.1016/j.copbio.2009.10.008.

13 Ghaderi D, et al. Production Platforms for Biotherapeutic Glycoproteins: Occurrence, Impact, and Challenges of Nonhuman Sialylation. *Biotechnol. Bioeng. Rev.* 28(1) 2012: 147–176.

14 Hossler P, Khattak SF, Jian Li Z. Optimal and Consistent Protein Glycosylation in Mammalian Cell Culture. *Glycobiology* 19(9) 2009: 936–949; doi:10.1093/glycob/cwp079.

15 Aricescu AR, Lu W, Jones EY. A Time and Cost-Efficient System for High-Level Protein Production in Mammalian Cells. *Acta Cryst. D62* 2006: 1243–1250; doi:10.1107/S0907444906029799.

16 Dalton A, Barton WA. Overexpression of Secreted Proteins from Mammalian Cells. *Protein Sci.* 23(5) 2014: 517–525; doi:10.1002/pro.2439.

17 Strasser R, Altmann F, Steinkellner H.

- Controlled Glycosylation of Plant-Produced Recombinant Proteins. *Curr. Opin. Biotechnol.* 30, 2014: 95–100; doi:10.1016/j.copbio.2014.06.008.
- 18 Steger K, et al. *Literature Review: CHO Versus HEK Cell Glycosylation*. Maxcyte: Gaithersburg, MD.
- 19 van Berkel PHC, et al. N-Linked Glycosylation Is an Important Parameter for Optimal Selection of Cell Lines Producing Biopharmaceutical Human IgG. *Biotechnol. Prog.* 25(1) 2009: 244–251; doi:10.1002/btpr.92.
- 20 Sheeley DM, Merrill DM, Taylor LCE. Characterization of Monoclonal Antibody Glycosylation: Comparison of Expression Systems and Identification of Terminal  $\alpha$ -Linked Galactose. *Anal. Biochem.* 247(1) 1997: 102–110.
- 21 Raju TS, Jordan RE. Galactosylation Variations in Marketed Therapeutic Antibodies. *MAbs* 4(3) 2012: 385–391.
- 22 Li F, et al. Cell Culture Processes for Monoclonal Antibody Production. *MAbs* 2(5) 2010: 466–477; doi:10.4161/mabs.2.5.12720.
- 23 Liu B, et al. The Availability of Glucose to CHO Cells Affects the Intracellular Lipid-Linked Oligosaccharide Distribution, Site Occupancy and the N-Glycosylation Profile of a Monoclonal Antibody. *J. Biotechnol.* 170, 2014: 17–27.
- 24 Surve T, Gadgil M. Manganese Increases High Mannose Glycoform on Monoclonal Antibody Expressed in CHO When Glucose Is Absent or Limiting: Implications for Alternate Sugars. *Biotechnol. Progress* 31(2) 2015: 460–467; doi:10.1002/btpr.2029.
- 25 Wong NS, et al. An Investigation of Intracellular Glycosylation Activities in CHO Cells: Effects of Nucleotide Sugar Precursor Feeding. *Biotechnol. Bioeng.* 107(2) 2010: 321–336; doi:10.1002/bit.22812.
- 26 Chen P, Harcum SW. Effects of Amino Acid Additions on Ammonium-Stressed CHO Cells. *J. Biotechnol.* 117(3) 2005: 277–286.
- 27 Borys MC, Linzer DI, Papoutsakis ET. Ammonia Affects the Glycosylation Patterns of Recombinant Mouse Placental Lactogen-I By Chinese Hamster Ovary Cells in a pH-Dependent Manner. *Biotechnol. Bioeng.* 43(6) 1994: 505–514; doi: 10.1002/bit.260430611.
- 28 McAfee A, Templeton N, Young J. Role of Chinese Hamster Ovary Central Carbon Metabolism in Controlling the Quality of Secreted Biotherapeutic Proteins. *Pharm. Bioprocess.* 2(1) 2014: 63–74; doi:10.4155/bpp.13.65.
- 29 Chen P, Harcum S. Effects of Elevated Ammonium on Glycosylation Gene Expression in CHO Cells. *Metabol. Eng.* 8(2) 2006: 123–132; doi:10.1016/j.jymben.2005.10.002.
- 30 Hong JK, Cho SM, Yoon SK. Substitution of Glutamine By Glutamate Enhances Production and Galactosylation of Recombinant IgG in Chinese Hamster Ovary Cells. *Appl. Microbiol. Biotechnol.* 88(4) 2010: 869–876.
- 31 Roth JM. Recombinant Tissue Plasminogen Activator for the Treatment of Acute Ischemic Stroke. *Proc. (Bayl. Univ. Med. Cent.)* 24(3) 2011: 257–259.
- 32 McCracken NA, Kowle R, Ouyang A. Control of Galactosylated Glycoforms Distribution in Cell Culture System. *Biotechnol. Prog.* 30(3) 2014: 547–553; doi:10.1002/btpr.1906.
- 33 Clincke M, et al. Effect of Iron Sources on the Glycosylation Macroheterogeneity of Human Recombinant IFN- $\gamma$  Produced By CHO Cells During Batch Process. *BMC Proceedings* 5, 2011: 114; doi:10.1186/1753-6561-5-S8-P114.
- 34 Hossler P, Racicot C, McDermott S. Targeted Shifting of Protein Glycosylation Profiles in Mammalian Cell Culture Through Media Supplementation of Cobalt. *J. Glycobiol.* 3, 2014: 1; doi:10.4172/2168-958X.1000108.
- 35 Yuk IH, et al. Effects of Copper on CHO Cells: Cellular Requirements and Product Quality Considerations. *Biotechnol. Prog.* 31(1) 2015: 226–238; doi:10.1002/btpr.2004.
- 36 Crowell CK, Grampp GE. *Production of Glycoproteins Using Manganese*. US Patent 7972810 B2, 2011.
- 37 Gramer MJ, et al. Modulation of Antibody Galactosylation Through Feeding of Uridine, Manganese Chloride, and Galactose. *Biotechnol. Bioeng.* 108(7) 2011: 1591–1602; doi: 10.1002/bit.23075.
- 38 Grainger RK, James DC. CHO Cell Line Specific Prediction and Control of Recombinant Monoclonal Antibody N-Glycosylation. *Biotechnol. Bioeng.* 110(11) 2013: 2970–2983; doi:10.1002/bit.24959.
- 39 DePalma A. Glycosylation: The Key to Quality. *Genet. Eng. Biotech. N.* 35(6) 2015.
- 40 Wilson S, et al. *Raw Material Characterization: Effect of Trace Metal Variability on Protein Glycosylation*. Sigma Aldrich Technical Poster. 

**Lori Fortin**, senior manager, R&D  
**Jonathan Fura**, manager, R&D  
**Hong Li**, formerly entrepreneurial scientist  
**Brian Beyer**, formerly director of bioprocess, R&D  
**Nandu Deorkar** is vice president of R&D at Avantor Inc., Bridgewater, NJ; nandu.deorkar@avantorsciences.com.

Reprinted with permission to Avantor from *BioProcess International*  
 January-February © 2018 Informa.